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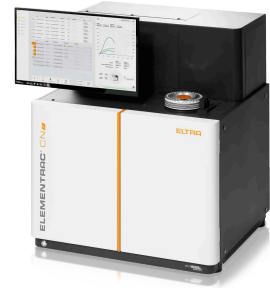
ELEMENTRAC CN-r

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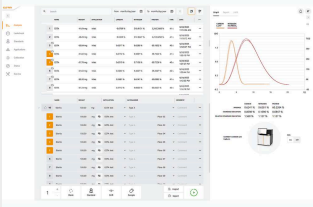
Step 1: Registration of the Sample in the Elements 4.0 Software

The sample is registered in the ELEMENTS 4.0 software. The weight is automatically transferred from the balance (step 1) – fast, safe, and error-free.

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The samples are conveniently placed into the carousel. The CN-r automatically detects the carousel and clearly assigns all samples – ensuring a smooth, safe, and automated process.

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
The ELEMENTRAC CN-r can be flexibly configured:


- | Nitrogen/Protein determination
- | Nitrogen/Protein and Carbon determination

While carbon is detected as CO₂, nitrogen determination is performed in elemental form using a highly sensitive thermal conductivity cell (TCD). Helium is used as the standard carrier gas (argon optional). This ensures maximum sensitivity in nitrogen analysis.

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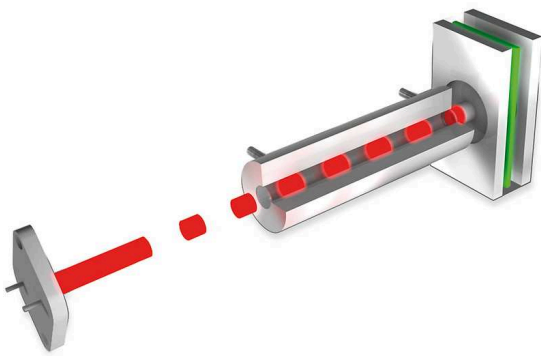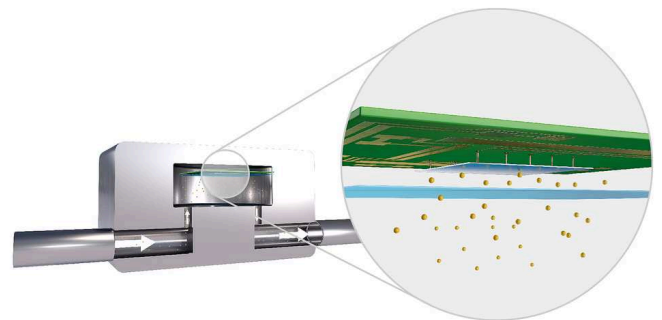All chemicals required for operation are ergonomically arranged at the front and safely hidden behind a door during routine use. Maintenance is especially user-friendly: combustion and reduction tubes can be replaced without tools in just a few minutes. This well-thought-out design reduces time expenditure, increases reliability, and ensures maximum reproducibility of your analyses.

TOUCHSCREEN & INTEGRATED PC

The ELEMENTRAC CN-r is equipped with a modern touchscreen that can be placed flexibly either directly on the device or conveniently on the desk. Thanks to the integrated PC, no additional hardware is required. The ELEMENTS 4.0 software can be intuitively operated via touchscreen – or classically with a keyboard and mouse. With the available USB ports, you can easily connect peripheral devices such as a keyboard, mouse, or a balance.

NETWORKING & DATA ACCESS

The CN-r can be easily integrated into your network. This allows you to transfer results directly into your LIMS or access the device software from your office PC. Whether recalculations, evaluations, or entering new samples: you always have full access and can complete your tasks easily and efficiently – regardless of the device's location.

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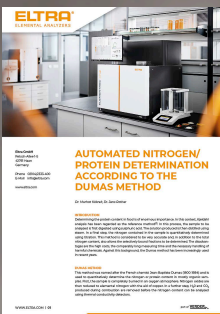
- | Determination of protein content for nutritional labeling and quality control
- | Evaluation of rice by-products (bran, husks) for animal feed regarding nutrient composition
- | Supporting breeding and research for improved rice varieties (higher protein content, better nutritional properties)

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ORDER DATA

ELEMENTRAC[®] CN-R

Measuring range for 300 mg sample weight

88200-4000		CN-r	0.007 – 58.3 % C (0,02 – 175 mg C abs.) 0.01 – 100 % N (0,03 – 300 mg N abs.)
88200-4001		CN-r	0.01 – 100 % N (0,03 – 300 mg N abs.)

REQUIRED ACCESSORIES

Linux based PC is already integrated, as well as 21,5 " Touchscreen monitor

88400-0645	Balance (resolution 0.0001 g)
------------	-------------------------------

REQUIRED CONSUMABLES / CHEMICALS FOR FIRST OPERATIONS

88500-0021	CN-r / N-r Starting kit
------------	-------------------------

90200		Anhydron
-------	-------------------------------------------------------------------------------------	----------

90210		Sodium hydroxide
-------	-------------------------------------------------------------------------------------	------------------

FURTHER OPTIONS AND CONSUMABLES

88400-0474	Tin foils, 36x36 mm, -HR-(200 pieces / packing unit)
------------	------------------------------------------------------

88400-0640	Copper Turnings, 200 g
------------	------------------------

88400-0638	N Catalyst, 50 g
------------	------------------

88400-0637	Large Ceramic Ash Crucibles-HR-(10 pieces / packing unit)
------------	-----------------------------------------------------------

88400-0636	Copper sticks-HR-(100 g/packing unit)
------------	---------------------------------------

88400-0635	Aluminium oxide pellets (Furnace reagent); 200 g
------------	--------------------------------------------------

88400-0508		Steel wool, 454 g
------------	-------------------------------------------------------------------------------------	-------------------

90330		Quartz wool (50 g)
-------	-------------------------------------------------------------------------------------	--------------------

90332  Glas wool (50 g)

88400-0644 Tinfoil, 51x51 mm (200 pieces / packing unit)

FURTHER HARDWARE TOOLS (SPATULA, TONGUES)

88400-0229  Tweezers (160 mm), curved, 1 piece

88400-0472  Tweezers (145 mm), straight, 1 piece

88400-0475  Set with 6 spatula and 1 tweezers for multiple weighing procedures

TYPICAL SPARE PARTS

88400-0006  Reagent glass tube 20x280 mm

11064-3002 Reagent tube 305x20

40000-3512 Combustion tube

CALIBRATION MATERIALS

PRIMARY SUBSTANCES

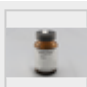
90710-3010  EDTA, 50 g

90710-3030 Succrose

90810  Calcium carbonate

COAL PREMIUM

92550-3010 Specified for C;H;N;S ; TGA

92550-3020  Specified for C;H;N;S ; TGA

92550-3040

Specified for C;H;N;S ; TGA

92550-3060

Specified for C;H;N;S ; TGA

COKE PREMIUM

92560-3010

Specified for C;H;N;S ; TGA

PET COKE PREMIUM

92570-3020

Specified for C;H;N;S ; TGA

92570-3020

Specified for C;H;N;S ; TGA