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CHS-580A

**ELTRA's CHS-580A is the ideal elemental analyzer for simultaneous determination of carbon, hydrogen and sulfur in organic samples. It features a resistance furnace with vertical ceramic tube and uses ceramic crucibles as sample carriers.**

An automatic sample loader allows for analysis of large numbers of samples without operator intervention. The sample loader is available in two versions to accommodate 36 or 130 crucibles. For reliable analysis of acidified samples (TOC analysis), both versions are available in an acid-resistance configuration.

The temperature can be set in steps of 1 °C to a maximum of 1,550 °C. Thanks to sample weights of up to 500 mg and more even inhomogeneous materials can be reliably analyzed. The analyzer can be equipped with up to four independent infrared cells for measuring carbon and sulfur concentrations, allowing for a great variety of applications.

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- | Fast & simultaneous elemental analysis of C, H, S, TOC and TIC
- | Configurable for single element (C or S) or multi-element (C/S or C/H/S) analysis
- | Wide concentration range from 0.005 – 100 %
- | High weights possible for elemental analysis of heterogeneous samples
- | Optional autoloader with 36 or 130 positions (acid-resistant version available)

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## OPERATION AND ANALYSIS PROCESS

Operation of the CHS-580 A elemental analyzer is simple, fast and convenient. The samples are weighed and registered in the software as manual analysis or analysis via autoloader. Typical sample weights are between 50 and 500 mg. After starting the analysis, the sample is introduced into the furnace and the combustion gases CO<sub>2</sub> and SO<sub>2</sub> are measured in up to four infrared measuring cells. The results of the elemental analysis can be exported via LIMS, report or text file.

*Logging the sample into the software and weighing*

*Placing sample in autoloader and analysis*

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### **CONFIGURATIONS AND OPTIONS**

ELTRA CHS-580 A analyzers can be configured to meet a wide range of customer requirements. C, H, S and C/H/S configurations are available with up to 2 infrared cells for each element to cover a wide working range. Robust measuring cells for the reliable determination of sulfur in very high concentrations are also available.

The CHS-580 A series is perfectly equipped for TOC analysis according to DIN EN 15936 with optional acid-resistant autosamplers, as well as the optional TIC module.

An optional autoloader with 130 crucible positions is available for analyses with high sample throughput.

*Analyzer with optional TIC module*

*Autoloader with 130 crucible positions*

*Autoloader with acid-resistant plastic crucible holders*

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In the CHS-580A analyzers the sample is burnt in an oxygen atmosphere at temperatures of up to 1,550 °C. The furnace temperature can be freely selected in steps of 1 °C. The combustion gasses (CO<sub>2</sub>, H<sub>2</sub>O, SO<sub>2</sub>) coming from the furnace pass through a dust filter. After the water vapor is chemically absorbed, the dried CO<sub>2</sub> and SO<sub>2</sub> gas is detected in the additional infrared cells. Depending on the configuration, it is possible to combine up to four infrared cells with different sensitivities.

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#### REQUIRED CONSUMABLES / CHEMICALS FOR FIRST OPERATIONS

88500-0006	Starter-kit for 500 analyses CHS 580A (500 crucibles, 50 g glass wool, 100 g Combsolid)
------------	---

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### CHEMICALS (FILLINGS FOR GLASS AND QUARTZ TUBES)

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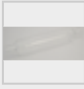
## TOOLS FOR OPERATION: SPATULAS, TWEEZERS, TONGS AND OTHER

88400-0229		Пинцет (160 мм), скругленный, 1 штука
88400-0472		Пинцет (145 мм), прямой, 1 штука
88400-0475		Set with 6 spatula and 1 tweezers
88400-0477		Weighing boat, 1 piece <span style="float: right;">for weighing and usage of granulates</span>
90145		Tongs for ceramic crucibles and boats, 220 mm, 1 piece
90146		Steel pot for preheated crucibles and used crucibles

## TOOLS FOR MAINTENANCE

35400-8000	Maintenance kit CS-580A
------------	-------------------------

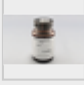
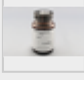
71010		Brush, 16 mm, 1 piece, for cleaning balance from dust
88400-0473		Powder funnel (plastics), 1 piece, for easy filling of chemical tubes
88400-0490		Rubber plug 29 x 35 x 30 mm, 1 piece, for sealing big glass tubes like 09090
88600-0026		Anhydrone filter tube

## CALIBRATION MATERIALS

**Calibration materials may show slight variations depending on the current lot.**

**To see the current certification please visit [www.ELTRA.com](http://www.ELTRA.com).**

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### COAL, PREMIUM, C/H/N/S, ASH, VOLATILE CONTENT, CERTIFIED

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**COKE, PREMIUM, C/H/N/S, ASH, VOLATILE CONTENT, CERTIFIED**

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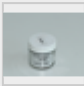
**PET COKE, PREMIUM, C/H/N/S, ASH, VOLATILE CONTENT CERTIFIED**

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90817-3003 Soil, 25 g, > 2 % C, S

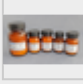
90817-3004 Soil, 25 g, > 2 % C, S

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PRIMARY SUBSTANCES FOR CALIBRATION

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LIQUID CALIBRATION MATERIAL

**Please note: Every analyzer requires PC, monitor, balance and some consumables (crucibles, chemicals) which have to be ordered separately**