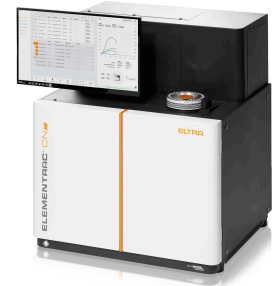
ANALIZOR DE CARBON / AZOT / PROTEINE

ELEMENTRAC CN-r

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

ENTRAC-CN

ANALIZOR DE CARBON / AZOT / PROTEINE
ELEMENTRAC CN-r

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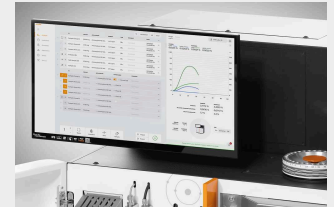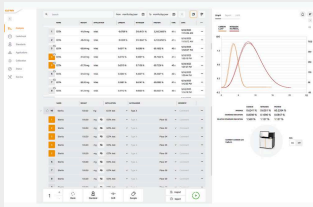Step 1: Registration of the Sample in the Elements 4.0 Software

The sample is registered in the ELEMENTS 4.0 software. The weight is automatically transferred from the balance (step 1) – fast, safe, and error-free.

Step 2: Weighing and Application

Whether a few milligrams or up to 1 gram: the ELEMENTRAC CN-r offers the highest precision with various sample types – solid or liquid. These are simply weighed into tin foil or tin capsules and reliably analyzed.

Pasul 3: Analiza

The samples are conveniently placed into the carousel. The CN-r automatically detects the carousel and clearly assigns all samples – ensuring a smooth, safe, and automated process.

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ANALIZOR DE CARBON / AZOT / PROTEINE
ELEMENTRAC CN-r

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
The ELEMENTRAC CN-r can be flexibly configured:


- | Nitrogen/Protein determination
- | Nitrogen/Protein and Carbon determination

While carbon is detected as CO₂, nitrogen determination is performed in elemental form using a highly sensitive thermal conductivity cell (TCD). Helium is used as the standard carrier gas (argon optional). This ensures maximum sensitivity in nitrogen analysis.

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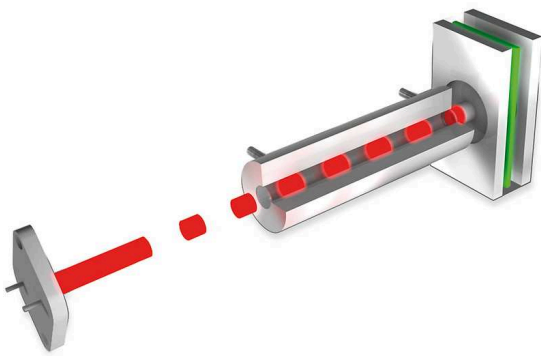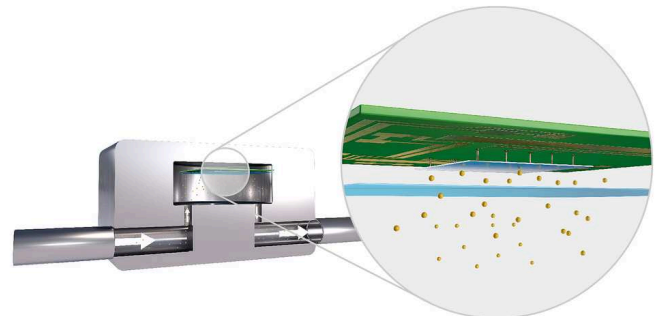ANALIZOR DE CARBON / AZOT / PROTEINE
ELEMENTRAC CN-r

SOLUȚII STANDARD INTEGRATE

All chemicals required for operation are ergonomically arranged at the front and safely hidden behind a door during routine use. Maintenance is especially user-friendly: combustion and reduction tubes can be replaced without tools in just a few minutes. This well-thought-out design reduces time expenditure, increases reliability, and ensures maximum reproducibility of your analyses.

TOUCHSCREEN & INTEGRATED PC

The ELEMENTRAC CN-r is equipped with a modern touchscreen that can be placed flexibly either directly on the device or conveniently on the desk. Thanks to the integrated PC, no additional hardware is required. The ELEMENTS 4.0 software can be intuitively operated via touchscreen – or classically with a keyboard and mouse. With the available USB ports, you can easily connect peripheral devices such as a keyboard, mouse, or a balance.

NETWORKING & DATA ACCESS

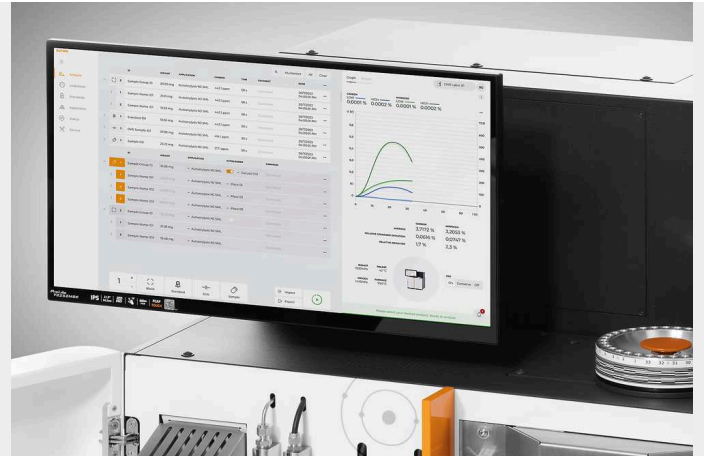
The CN-r can be easily integrated into your network. This allows you to transfer results directly into your LIMS or access the device software from your office PC. Whether recalculations, evaluations, or entering new samples: you always have full access and can complete your tasks easily and efficiently – regardless of the device's location.

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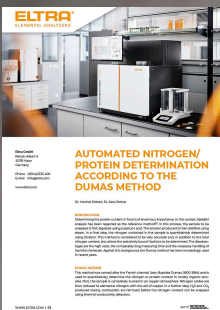
ANALIZOR DE CARBON / AZOT / PROTEINE
ELEMENTRAC CN-r

SOFTWARE

The comprehensive ELEMENTS software is an essential component of all elemental analyzers of the ELEMENTRAC generation.

A central window (analysis and results) is the starting point from which all functionalities required for the daily routine are easily accessible. From here it is possible to group and export analyzed samples or to register and analyze new samples. The user can access various subordinate functionalities such as application settings, calibration, diagnostics, audit trail, user group management or status and is FDA 21 CFR Part 11 compliant.

In addition, thanks to the network integrity and web browser architecture, you can access the device from any computer in the company network and have all data available at any time.

APPLICATION NOTES

PROTEIN DETERMINATION IN MILK WITH DUMAS METHOD

The Dumas method is a valid alternative for determining protein content in milk, providing quick and accurate results without the use of harmful chemicals. Its efficiency and safety

make it increasingly popular compared to the Kjeldahl method. Read our application note...

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
RICE

- | Determination of protein content for nutritional labeling and quality control
- | Evaluation of rice by-products (bran, husks) for animal feed regarding nutrient composition
- | Supporting breeding and research for improved rice varieties (higher protein content, better nutritional properties)

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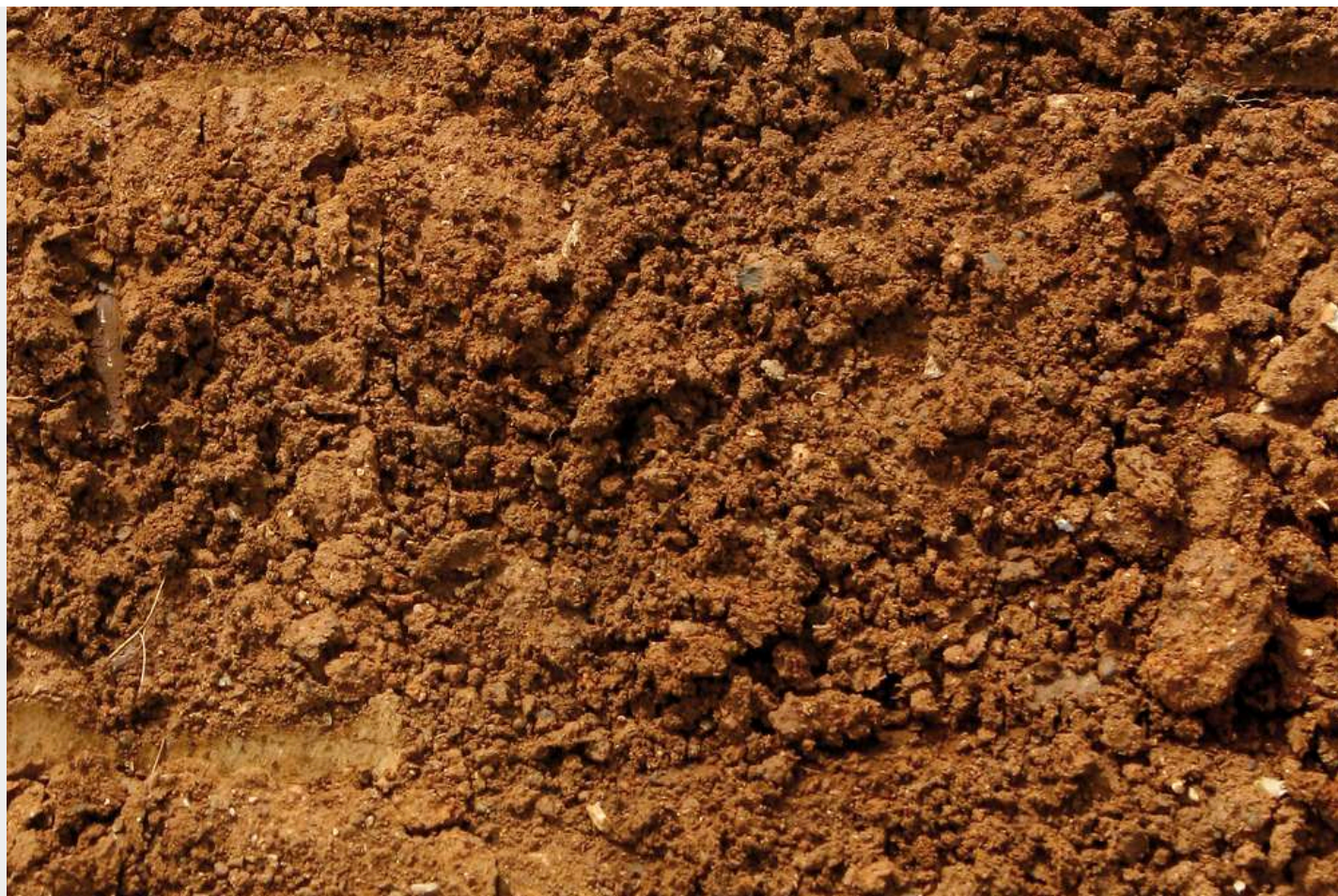MILK & DAIRY PRODUCTS

- | Analysis of solid or liquid samples such as cheese, blue cheese, yogurt, and milk
- | Determination of protein content for nutritional labeling and quality control
- | The ELEMENTRAC CN-r meets or exceeds the requirements of all common international standards, such as DIN EN ISO 14891

CARNE & PRODUSE DIN CARNE

- | Determination of protein content for nutritional labeling and quality control
- | Verification of meat authenticity and prevention of food fraud
- | Monitoring protein levels in processed meat products

COAL

- | Measurement of nitrogen content to assess quality
- | Control of nitrogen levels in coke production
- | Optimization of combustion processes and reduction of emissions

FERTILIZATORI

- | Determination of nitrogen concentration to ensure fertilizer efficiency
- | Quality control and compliance with agricultural standards
- | Evaluation of nutrient availability for improved crop yield

GRÂNE & SEMINȚE DE ULEI

- | Protein determination for nutritional evaluation and feed production
- | Quality assessment in oilseed processing (soy, sunflower, rapeseed, etc.)
- | Supporting research on oilseed breeding and agricultural development

ANALIZOR DE CARBON / AZOT / PROTEINE
ELEMENTRAC CN-r

DATE TEHNICE

Measurement ranges	Nitrogen: 0.03 mg to 300 mg Carbon: 0.02 mg to 175 mg
Typical weights	Up to 1.0 g, 0.5 g nominal
Analysis time	< 3 min
Detectors	Thermal conductivity detector (N) NDIR (C)
Typical samples	Food, feed stuff, fuels (coal and coke), environmental samples and fertilizers
Reagents	Magnesium perchlorate, sodium hydroxide, copper, aluminum oxide
Sursă de alimentare	230 VAC ± 10 %, 50/60 Hz
Gaze necesare	Oxygen (99.999% purity) Helium (99.996% purity) Compressed air (oil and water free)
Configurations	CN-r, N-r
Greutate	135 kg
Dimensiuni (W x H x D)	(W x H x D) 75,7 x 78,5 x 62 cm

www.eltra.com/cn-r

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ELEMENTRAC[®] CN-R

Measuring range for 300 mg sample weight

88200-4000		CN-r	0.007 – 58.3 % C (0,02 – 175 mg C abs.) 0.01 – 100 % N (0,03 – 300 mg N abs.)
88200-4001		CN-r	0.01 – 100 % N (0,03 – 300 mg N abs.)

REQUIRED ACCESSORIES

Linux based PC is already integrated, as well as 21,5 " Touchscreen monitor

88400-0645	Balance (resolution 0.0001 g)
------------	-------------------------------

REQUIRED CONSUMABLES / CHEMICALS FOR FIRST OPERATIONS

88500-0021	CN-r / N-r Starting kit
------------	-------------------------

90200		Anhydron
-------	---	----------

90210		Sodium hydroxide
-------	---	------------------

FURTHER OPTIONS AND CONSUMABLES

88400-0474	Tin foils, 36x36 mm, -HR-(200 pieces / packing unit)
------------	--

88400-0640	Copper Turnings, 200 g
------------	------------------------

88400-0638	N Catalyst, 50 g
------------	------------------

88400-0637	Large Ceramic Ash Crucibles-HR-(10 pieces / packing unit)
------------	---

88400-0636	Copper sticks-HR-(100 g/packing unit)
------------	---------------------------------------

88400-0635	Aluminium oxide pellets (Furnace reagent); 200 g
------------	--

88400-0508		Steel wool, 454 g
------------	---	-------------------

90330		Quartz wool (50 g)
-------	---	--------------------

90332

Glas wool (50 g)

88400-0644

Tinfoil, 51x51 mm (200 pieces / packing unit)

FURTHER HARDWARE TOOLS (SPATULA, TONGUES)

88400-0229

Tweezers (160 mm), curved, 1 piece

88400-0472

Tweezers (145 mm), straight, 1 piece

88400-0475

Set with 6 spatula and 1 tweezers for multiple weighing procedures

TYPICAL SPARE PARTS

88400-0006

Reagent glass tube 20x280 mm

11064-3002

Reagent tube 305x20

40000-3512

Combustion tube

CALIBRATION MATERIALS

PRIMARY SUBSTANCES

90710-3010

EDTA, 50 g

90710-3030

Succrose

90810

Calcium carbonate

COAL PREMIUM

92550-3010

Specified for C;H;N;S ; TGA

92550-3020

Specified for C;H;N;S ; TGA

92550-3040

Specified for C;H;N;S ; TGA

92550-3060

Specified for C;H;N;S ; TGA

COKE PREMIUM

92560-3010

Specified for C;H;N;S ; TGA

PET COKE PREMIUM

92570-3020

Specified for C;H;N;S ; TGA

92570-3020

Specified for C;H;N;S ; TGA